

FLIGHT ACT OFFICER LOG	DAY	REV	PG
SITE/ACQ/LOS	NOTES		
To - Biomed			
cc - Mike Whittle, Ray McKinney			
A clean up of M074 is needed - The following actions should be taken -			
1. Whatever is required to obtain Cal data performed some 4 days ago -	AETN/EVC		
2. Changeout of WMC SMMD electronic package - Suggest the following procedure:			
a) remove + replace zero-G connector (pwr off)			
b) turn pwr. switch to mass + temp. noting presence or absence of power supply squeal + lights (power on)			
c) Assuming unit has failed remove W.M.C. electronics and replace with W.R. electronics			
[changeout procedures must be generated]			
d) perform calibration per C/L			
W. Thornton			
Note - no fecal or mass data until all these procedures have been performed -			